

Cleaning the Tissue Culture Hood

Note: The tissue culture hood should only contain items that are being worked with - the pipettor, suction tube etc. All plastics and other equipment should be kept outside the hood. The hood is not a storage area!

Once a week:

1. Wear nitrile gloves, lab-coat and goggles.
2. Take all items out of the hood.
3. With a solution that is 1-part Spitfire-5 parts water, wash the surface and walls of the hood.
4. Let washed areas dry. Then rinse with a clean sponge and water.
5. Put items back in hood when finished.

Once a month:

1. Wear nitrile gloves, lab-coat and goggles.
2. Take all the items out of the hood.
3. Open the hood door and take out the hood benchtop.
4. Wash and rinse the inside of the hood with Spitfire solution as above